



Syrris Ltd.
27 Jarman Way, Royston, Hertfordshire
SG8 5HW, United Kingdom
T: +44 (0)1763 242555
F: +44 (0)1763 242992
E: info@syrris.com
W: www.syrris.com

Syrris Limited
Registered office as address
Company No. 4257809

Subject	Circulator Cooling Water Specification
Date	17-NOV-2011
Author	SV

Please find below the specification of the water used in water cooled circulator units.

Failure to ensure the water specification is met will result in poor performance from the circulator. In a worst case scenario the circulator will simply show an error message and not work.

	Huber	Julabo
Cooling Water Temperature	15DegC	<20DegC
Maximum Input Pressure	6 bar	6 bar
Pressure Difference (In-Out)	3 bar +	3.5 to 6 bar
pH at 25DegC	7	7 to 8.5
Suspended Matter	-	<30mg/Litre
Size of suspended matter	-	<0.1mm

This information was collected from the Huber and Julabo service and support departments in Germany.

Please contact Support@Syrris.com if you require further information.